

General Cell Culture Protocol

Feeding Cells:

- 1) Clean and sterilize the hood using 70% ethanol
- 2) In water bath set for 37 °C, warm for 1 hour Complete Media and other necessary media
- 3) To Feed Cells:
 - a) Use pre-sterilized Louis Pasteur pipette attached to vacuum tubing to remove media from plate/flask.
 - b) Quickly add Complete Media to plate using sterile Falcon Pipet with Pipet-Aid, only feeding 1-3 plates of same cells at a time to avoid drying cells.
 - c) On the plate cover write the date fed and initial
 - d) Return plate to incubator